



**ken-a-vision®**  
KNOWLEDGE THROUGH VISION

# PupilCAM


## Instruction Manual

1401KRM

1401KEM

1402KRM

1402KEM

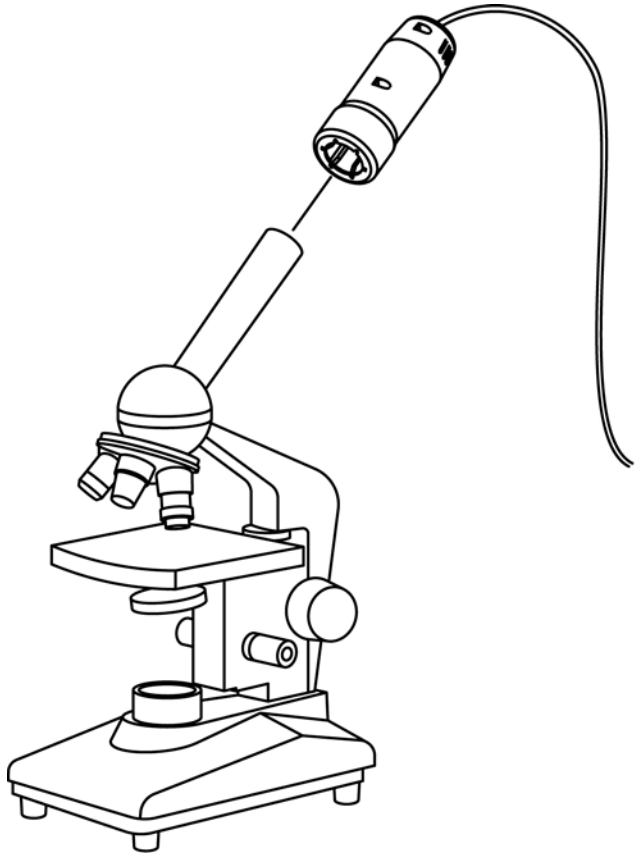

Thank you for your purchase of a Ken-A-Vision PupilCAM! This high-resolution camera captures microscopic images and displays them on your computer with full screen viewing. The pre-attached cable provides a direct interface to your equipment. It's ready to go, no assembly required!

## Connecting The USB Cable to Your Computer

Prior to connecting the USB cable, you must first install the supplied software.

Pre-attached cabling makes it easy to connect the USB cable to your computer. Simply insert the rectangular end of the USB cable to the USB port of your computer.

Note: The USB port will supply ample power to the the PupilCAM<sup>®</sup>. No power additional power supply is required.

## Connecting to a Microscope

The PupilCAM<sup>®</sup> is designed exclusively to work with a microscope (compound or stereoscope). A variable eyepiece adapter is built-in to the camera.

Gently twist the PupilCAM<sup>®</sup> as you push it down onto the microscope.

1401KRM/1402KRM models feature a rubber adapter that fits snugly over 28mm microscope eyepieces.

1401KEM/1402KEM models feature a built-in 10x eyepiece that is inserted into 23mm microscope eyepiece tubes or 30mm eyepiece tubes with the included adapter.

All focusing should be done using the microscope focus knobs.

To learn more about other Ken-A-Vision products  
and services please visit our website at  
[www.ken-a-vision.com](http://www.ken-a-vision.com)  
Or call  
1-816-353-4787

