


# Quick Start Guide:

Ken-A-Vision FlexCam® 2 Document Camera,  
910-171-200




## Step 1 Unpack Box

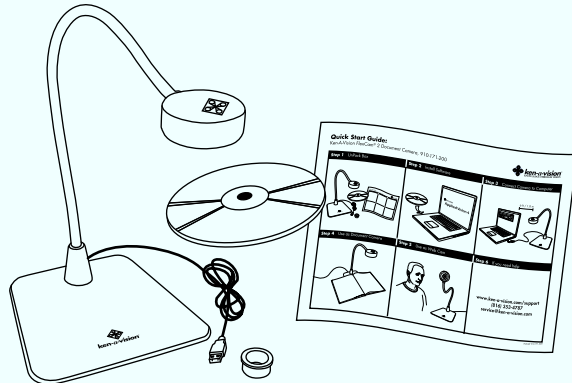


1. Camera  
(head, gooseneck body, base, eyepiece adapter, pre-attached USB cable)
2. CD with Applied Vision 4 (AV4) software  
(also downloadable at [www.ken-a-vision.com](http://www.ken-a-vision.com))
3. This Quick Start Guide


## Step 2 Install Software




1. Insert CD into your computer (PC or Mac).
2. Follow the onscreen prompts during installation.
3. Once installed, the AV4 icon  will appear on your desktop.

## Step 3 Connect Camera to Computer

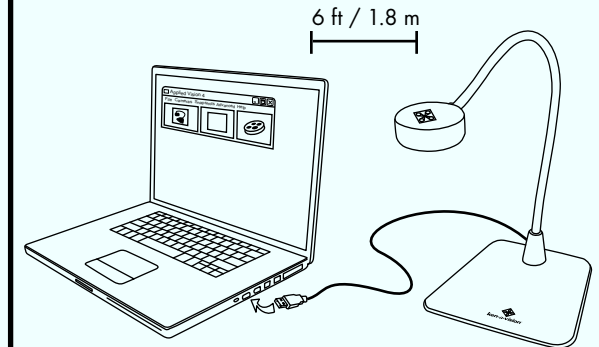


1. Position camera within 6 feet / 1.8 meters of the computer.
2. Connect the USB cable that is attached to the camera to your computer (PC or Mac).
3. Click on AV4 icon  to launch feature-rich applications.
4. The camera will launch any program that has a video interface (Skype, Windows Media Player).

## Step 4 Use as Document Camera

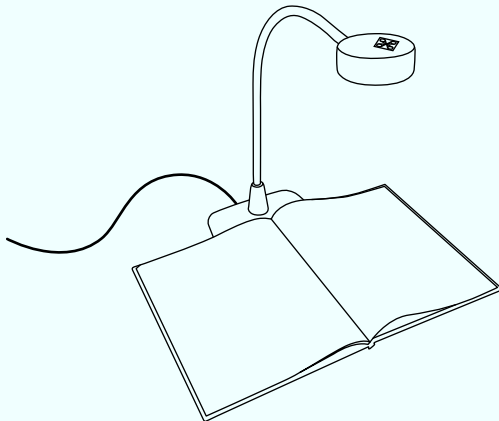


1. Place book, document or 3D object under the camera head. The base is flat and stable so you can place the item on top of it.
2. Move the camera head to adjust the image size and allow camera to focus.

## Step 5 Use as Web Cam

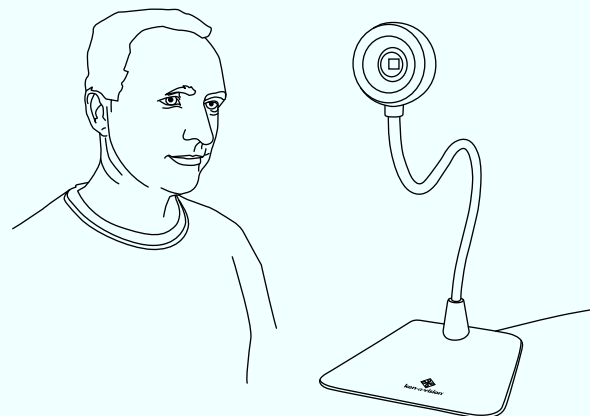

1. Flip up the camera head and aim at subject.
2. To rotate the image, use your keyboard and click on the rotate buttons.

## Step 6 If you need help

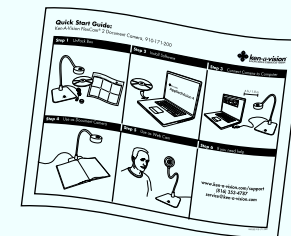

[www.ken-a-vision.com/support](http://www.ken-a-vision.com/support)  
**(816) 353-4787**  
[service@ken-a-vision.com](mailto:service@ken-a-vision.com)


1. **Visit:** [www.ken-a-vision.com](http://www.ken-a-vision.com) under Support
2. **Call:** (816) 353-4787
3. **Email:** [service@ken-a-vision.com](mailto:service@ken-a-vision.com)

# Quick Start Guide:

Ken-A-Vision FlexCam® 2 Document Camera,  
910-171-200




## Step 1 Unpack Box

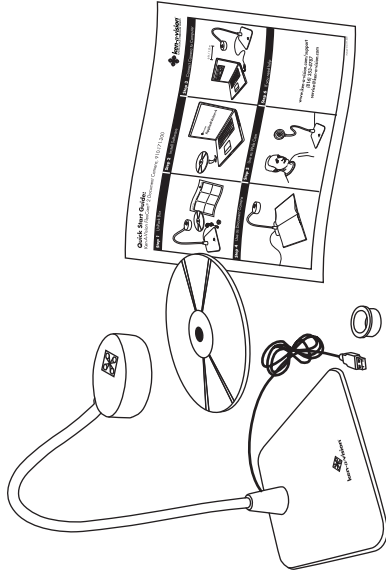


1. Camera (head, gooseneck body, base, eyepiece adapter, pre-attached USB cable)
2. CD with Applied Vision 4 (AV4) software (also downloadable at [www.ken-a-vision.com](http://www.ken-a-vision.com))
3. This Quick Start Guide

## Step 4 Use as Document Camera

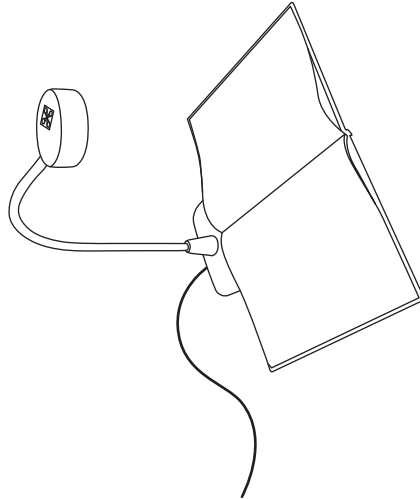


1. Place book, document or 3D object under the camera head. The base is flat and stable so you can place the item on top of it.
2. Move the camera head to adjust the image size and allow camera to focus.

## Step 2 Install Software

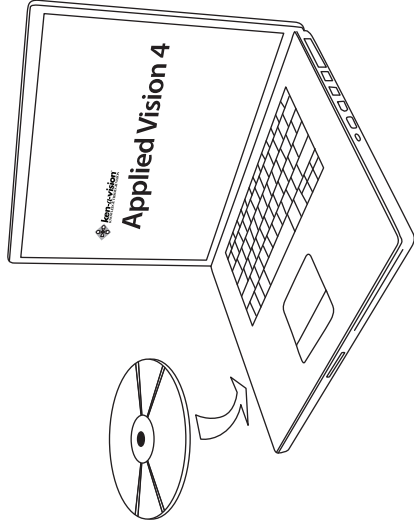


1. Insert CD into your computer (PC or Mac).
2. Follow the onscreen prompts during installation.
3. Once installed, the AV4 icon  will appear on your desktop.


## Step 5 Use as Web Cam

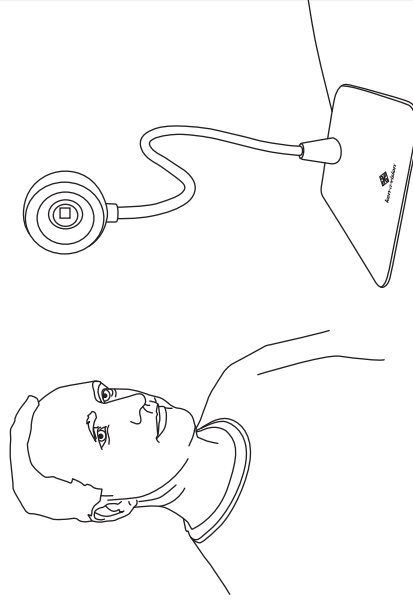


1. Flip up the camera head and aim at subject.
2. To rotate the image, use your keyboard and click on the rotate buttons.

## Step 3 Connect Camera to Computer

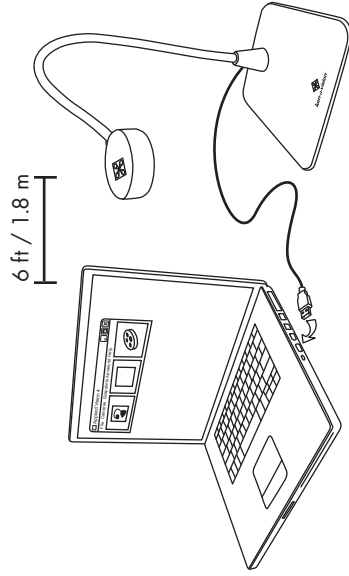

1. Position camera within 6 feet / 1.8 meters of the computer.
2. Connect the USB cable that is attached to the camera to your computer (PC or Mac).
3. Click on AV4 icon  to launch feature-rich applications.
4. The camera will launch any program that has a video interface (Skype, Windows Media Player).

## Step 6 If you need help

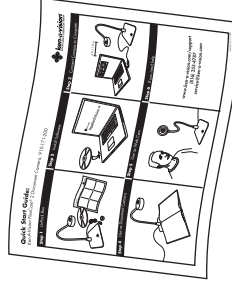

[www.ken-a-vision.com/support](http://www.ken-a-vision.com/support)  
**(816) 353-4787**  
**service@ken-a-vision.com**

1. **Visit:** [www.ken-a-vision.com](http://www.ken-a-vision.com) under Support
2. **Call:** (816) 353-4787
3. **Email:** [service@ken-a-vision.com](mailto:service@ken-a-vision.com)